AVANTOR VWR OX 4110 H OPERATING MANUAL

Available languages

-  English, Page 3
-  Deutsch, Seite 8
-  Español, Página 11
-  Français, Page 12
-  Italiano, Pagina 14
-  Português, Página 16
-  Čeština, Strana 18
-  Dansk, Side 20
-  ελληνικά, Σελίδα 22
-  Suomi, Sivu 24
-  Magyar, oldal 26
-  Dutch, Pagina 28
-  Norsk, Side 30
-  Polski, Strona 32
-  Română, Pagina 34
-  Русский, Страница 36
-  Slovenčina, Strana 38
-  Svenska, Sida 40
-  Türkçe, Sayfa 42

Quick Links

Deutsch

Table of Contents

[Table of Contents](#)

[English](#)

[Deutsch](#)

[Español](#)

[Français](#)

[Italiano](#)

[Português](#)

[CS Český](#)

[Dansk](#)

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[Svenska](#)

[TR Türkçe](#)

[Manual Versions in Other Languages](#)

[Other Manuals](#)[Lib Projects](#)

Operating manual

VWR® OX 4110 H - D.O. Meter

EU cat. no	NA cat. no
664-0176	-

CE

UK
CA

Legal Address of Manufacturer:

Europe

VWR International BV
Researchpark Haasrode 2020
Geldenaaksebaan 464
B-3001 Leuven
+32 16 385011
<http://be.vwr.com>

United States

VWR International LLC
100 Matsonford Rd
Radnor, PA 19087
+1 800-932-5000
www.vwr.com

Importer to UK:

VWR International Ltd
Hunter Boulevard, Magna Park
Lutterworth, Leicestershire, LE17 4XN
<http://uk.vwr.com>

Country of origin: Germany.

Contents

EN English	4
DE Deutsch.....	6
ES Español.....	8
FR Français	10
IT Italiano	12
PT Português	14
CS Český	16
DA Dansk	18
EL Ελληνικά	20
FI Suomi	22
HU Magyar	24
NL Nederlands	26
NO Norsk	28
PL Polski.....	30
RO Română	32
RU Русский	34
SK Slovenský	36
SV Svenska	38
TR Türkçe	40

OX 4110 H

Overview

The OX 4110 H meter enables you to perform measurements (D.O.) quickly and reliably.

The OX 4110 H meter provides a maximum degree of operating comfort, reliability and measuring certainty.

Keypad

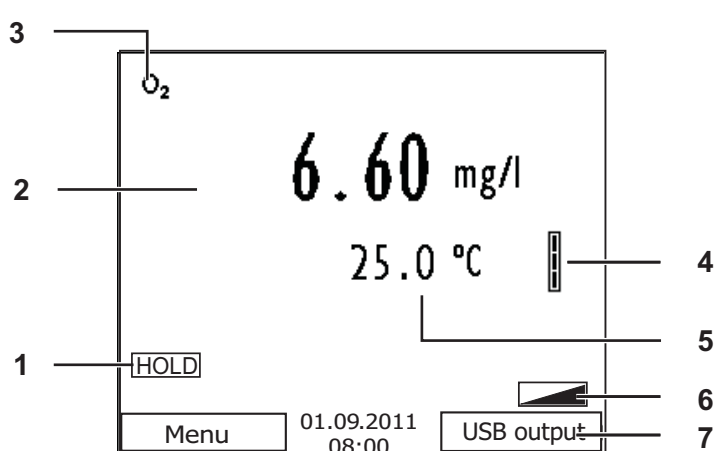
Keys are indicated by brackets <..>.

The key symbol (e.g. <OK>) generally indicates a short keystroke (under 2 sec).

A long keystroke (approx. 2 sec) is indicated by the underscore behind the key symbol (e.g. <OK__>).


<F1>:	Situation dependent function (is shown on the display) e.g.:
<F1__>:	<F1>/[Menu]: Opens the menu for measurement settings
<F2>:	<F1__>/[Menu]: Opens the menu for system settings
<F2__>:	
<On/Off>:	Switches the meter on or off
<MODE>:	Selects the measured parameter
<CAL>:	Calls up the calibration procedure
<CAL__>:	Displays the calibration data
<STR>:	Saves a measured value manually
<STR__> :	Opens the menu for the automatic save function
<RCL>:	Displays the manually stored measured values
<RCL__>:	Displays the automatically stored measured values
<▲><▼>:	Menu control, navigation
<OK>:	Opens the menu for measurement settings / confirms entries
<OK__>:	Opens the menu for system settings
<HOLD>:	Freezes the measured value (HOLD function)

Display

1	Status information
2	Measured value (with unit)
3	Measured parameter
4	Sensor symbol (calibration evaluation)
5	Measured temperature (with unit)
6	Further status information
7	Softkeys and date/time

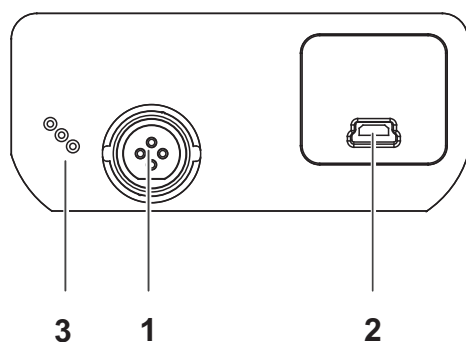Jack field /connections

1	D.O. sensor
2	USB B interface (device)
3	Service interface

Initial
commissioning

- 1 Insert the supplied batteries.
Make sure that the poles are in the correct position.
- 2 Switch on the meter with the <On/Off> key.

Caution

Follow the safety instructions of the used sensors.
The operating manuals of the sensors are available on the supplied CD.

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Local VWR offices in Europe and Asia Pacific

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India

VWR Lab Products Private Limited
No.139. BDA Industrial Suburb,
6th Main, Tumkur Road, Peenya Post,
Bangalore, India – 560058
Tel.: +91-80-28078400
Email: vwr_india@vwr.com

Ireland / Northern Ireland

VWR International Ltd / VWR
International (Northern Ireland) Ltd
Orion Business Campus
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: +353 01 88 22 222
Email sales.ie@vwr.com

Italy

VWR International S.r.l.
Via San Giusto 85
20153 Milano (MI)
Tel.: +39 02-3320311/02-487791
Email: info.it@vwr.com

Korea

VWR International ~
17, Daehak 4-ro, Yeongtong-gu,
Suwon-si, Gyeonggi-do
Tel.: +82 31 645 7256
saleskorea@avantorsciences.com

Middle East & Africa

VWR International FZ-LLC
DSP Laboratory Complex
125, Floor 01
Dubai, United Arab Emirates
Tel.: +971 4 5573271
info.mea@vwr.com

The Netherlands

VWR International B.V.
Postbus 8198
1005 AD Amsterdam
Tel.: +31 020 4808 400
Email: info.nl@vwr.com

Norway

VWR International AS
Brynsalleen 4
0667 Oslo
Tel.: +47 22 90 00 00
Email: info.no@vwr.com

Poland

VWR International Sp. z o.o.
Limbowa 5
80-175 Gdansk
Tel.: +48 058 32 38 210
Email: info.pl@vwr.com

Portugal

VWR International –
Material de Laboratório, Lda
Centro Empresarial de Alfragide
Rua da Indústria, nº 6
2610-088 Amadora
Tel.: +351 21 3600 770
Email: info.pt@vwr.com

Singapore

VWR Singapore Pte Ltd
18 Gul Drive
Singapore 629468
Tel: +65 6505 0760
Email: sales.sg@vwr.com

Spain

VWR International Eurolab S.L.
C/ Tecnología 5-17
A-7 Llinars Park
08450 - Llinars del Vallès
Barcelona
Tel.: +34 902 222 897
Email: info.es@vwr.com

Sweden

VWR International AB
Fagerstagatan 18a
163 94 Stockholm
Tel.: +46 08 621 34 00
Email: info.se@vwr.com

Switzerland

VWR International GmbH
Lerzenstrasse 16/18
8953 Dietikon
Tel.: +41 044 745 13 13
Email: info.ch@vwr.com

UK

VWR International Ltd
Customer Service Centre
Hunter Boulevard
Magna Park
Lutterworth
Leicestershire
LE17 4XN
Tel.: +44 (0) 800 22 33 44
Email: uksales@vwr.com

Manual Versions in Other Languages

Deutsch

Español

Mexicano

Français

Other ManualsLib Projects

www.manualslib.com

www.manualslib.de

www.manualslib.es

www.manualslib.fr

www.manualslib.nl

www.manualslib.mx

www.manualslib.tech 30+ Languages